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Liquids Analysis					
CFU/side	<25	25-50	50-200	>200	Not Evaluable
CFU/ml	10 ³	10 ⁴	10 ⁵	10 ⁶	10 ⁷

* 1 CFU/side = 1 / 1.25 x 10⁴ CFU/10m²

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MEDIUM		<i>Escherichia coli</i>	<i>Proteus spp</i>	<i>Klebsiella spp</i>	<i>Pseudomonas spp</i>	<i>Streptococcus spp</i>	<i>Staphylococcus spp</i>	Yeasts/Moulds
V.R.B.G. AGAR	Colonies appearance	RED-PURPLE	RED-PURPLE	RED-PURPLE	COLORLESS	INHIBITION	INHIBITION	INHIBITION
	Medium colour	RED-PURPLE	RED-PURPLE	RED-PURPLE	RED			
V.R.B.L. AGAR	Colonies appearance	RED-PURPLE	COLORLESS	RED-PURPLE	COLORLESS	INHIBITION	INHIBITION	
	Medium colour	RED-PURPLE	COLORLESS	RED-PURPLE	RED			
ROSE BENGAL CAF AGAR	Colonies appearance	INHIBITION	INHIBITION	INHIBITION	PARTIAL INHIBITION	INHIBITION	INHIBITION	PINK -WHITE
	Medium colour							PINK
VOGEL JOHNSON AGAR	Colonies appearance	INHIBITION	PARTIAL INHIBITION	INHIBITION	INHIBITION	INHIBITION	BLACK	
	Medium colour						YELLOW	
BILE ESCULIN AGAR	Colonies appearance	INHIBITION	PARTIAL INHIBITION	PARTIAL INHIBITION	INHIBITION	WHITE	INHIBITION	
	Medium colour					BLACK		
CETRIMIDE AGAR	Colonies appearance				GREEN			
	Medium colour				BIANCO			
PLATE COUNT AGAR + TTC	Colonies appearance	RED	RED	RED	RED	RED	RED	RED
	Medium colour	YELLOW	YELLOW	YELLOW	YELLOW	YELLOW	YELLOW	YELLOW

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