



MicroSnapTM | Enhanced Nutrient Broth

MicroSnap Enhanced Nutrient Broth is a ready-to-use, non-selective protein solution. The 9mL broth volume enables low level enumeration of product samples with MicroSnap test devices in 6 – 8 hours. The broth is specifically designed for 1mL milk samples and is compatible with MicroSnap Total, Coliform and *E. coli* test devices. MicroSnap Enhanced Nutrient Broth enables enumeration and low level detection (1CFU/mL milk) in 8 hours or less in a simple, low cost system.



Find support documents, instructional videos, and more at www.hygiena.com

Benefits:

- Low level enumeration in 6-8 hours (1 CFU/mL milk)
- Compatible with multiple detection tests (TVC, Coliform, and *E. coli*)
- 9 mL broth volume reduces interference from physical components in the sample
- Robust chemistry reduces interference from chemical interactions
- 12 month shelf life at refrigerated temperatures (2 – 8 °C)



Compatible Test Devices:

MicroSnap Total

With Enhanced Nutrient Broth, MicroSnap Total delivers TVC test results in 6 – 8 hours, giving you the advantage of knowing microbial levels in the same working day or shift. Presence/absence results are available within 8 hours. Enumeration results are available in as early as 6 hours.

MicroSnap Coliform & *E. coli*

MicroSnap Coliform and *E. coli* are rapid tests for detection and enumeration of Coliform and *E. coli* bacteria. With Enhanced Nutrient Broth, presence/absence results for milk samples are available within 8 hours.

Catalog No.	Description	Quantity
MS1-N-BROTH-9ML	MicroSnap Enhanced Nutrient Broth (9mL)	100
ENSURE	EnSURE Monitoring System with SureTrend Software	1
MS2-TOTAL	MicroSnap Total (TVC) Detection Device	100
MS2-ECOLI	MicroSnap <i>E. coli</i> Detection Device	100
MS2-COLIFORM	MicroSnap Coliform Detection Device	100



MicroSnapTM | Enhanced EB Broth

MicroSnap Enhanced EB Broth is a ready-to-use, non-selective protein solution. The 9mL broth volume enables low level enumeration of product samples with MicroSnap test devices in 6 – 8 hours. The broth is specifically designed for 1mL milk samples and is compatible with MicroSnap EB, Coliform, and *E. coli* test devices. MicroSnap Enhanced EB Broth enables enumeration and low level detection (1CFU/mL milk) in 8 hours or less in a simple, low cost system.



Find support documents, instructional videos, and more at www.hygiena.com

Benefits:

- Low level enumeration in 6-8 hours (1 CFU/mL milk)
- Compatible with multiple detection tests (EB, Coliform, and *E. coli*)
- 9 mL broth volume reduces interference from physical components in the sample
- Robust chemistry reduces interference from chemical interactions
- 12 month shelf life at refrigerated temperatures (2 – 8 °C)



Compatible Test Devices:

MicroSnap EB

With Enhanced EB Broth, MicroSnap EB delivers Enterobacteriaceae test results in 6 – 8 hours, giving you the advantage of knowing microbial levels in the same working day or shift. Presence/absence results are available within 8 hours. Enumeration results are available in as early as 6 hours.

MicroSnap Coliform & *E. coli*

MicroSnap Coliform and *E. coli* are rapid tests for detection and enumeration of Coliform and *E. coli* bacteria. With Enhanced EB Broth, presence/absence results for milk samples are available within 8 hours.

Catalog No.	Description	Quantity
MS1-EB-BROTH-9ML	MicroSnap Enhanced EB Broth (9mL)	100
ENSURE	EnSURE Monitoring System with SureTrend Software	1
MS2-EB	MicroSnap EB (Enterobacteriaceae) Detection Device	100
MS2-ECOLI	MicroSnap <i>E. coli</i> Detection Device	100
MS2-COLIFORM	MicroSnap Coliform Detection Device	100