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
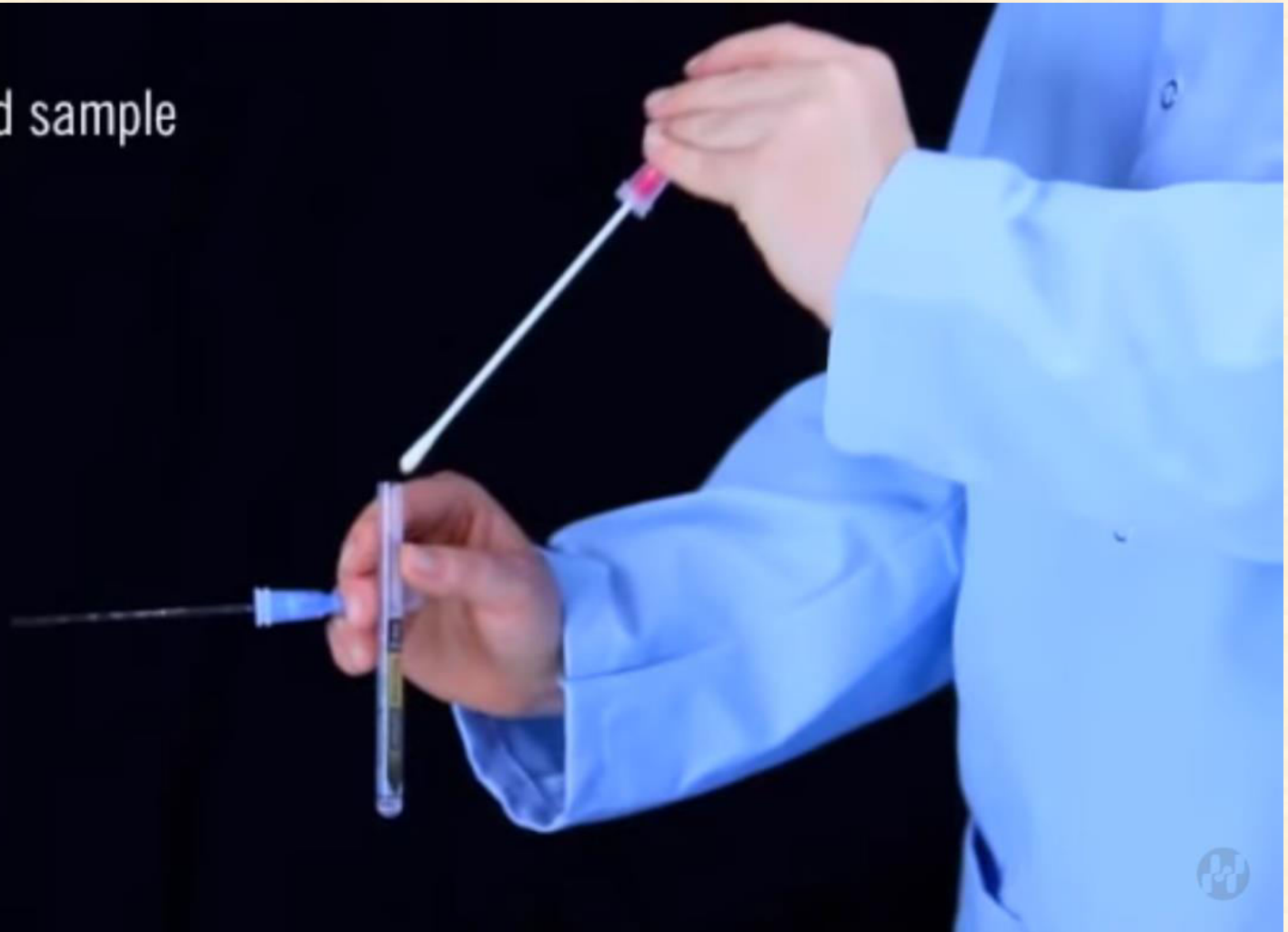
Pipette enriched sample
to fill line.

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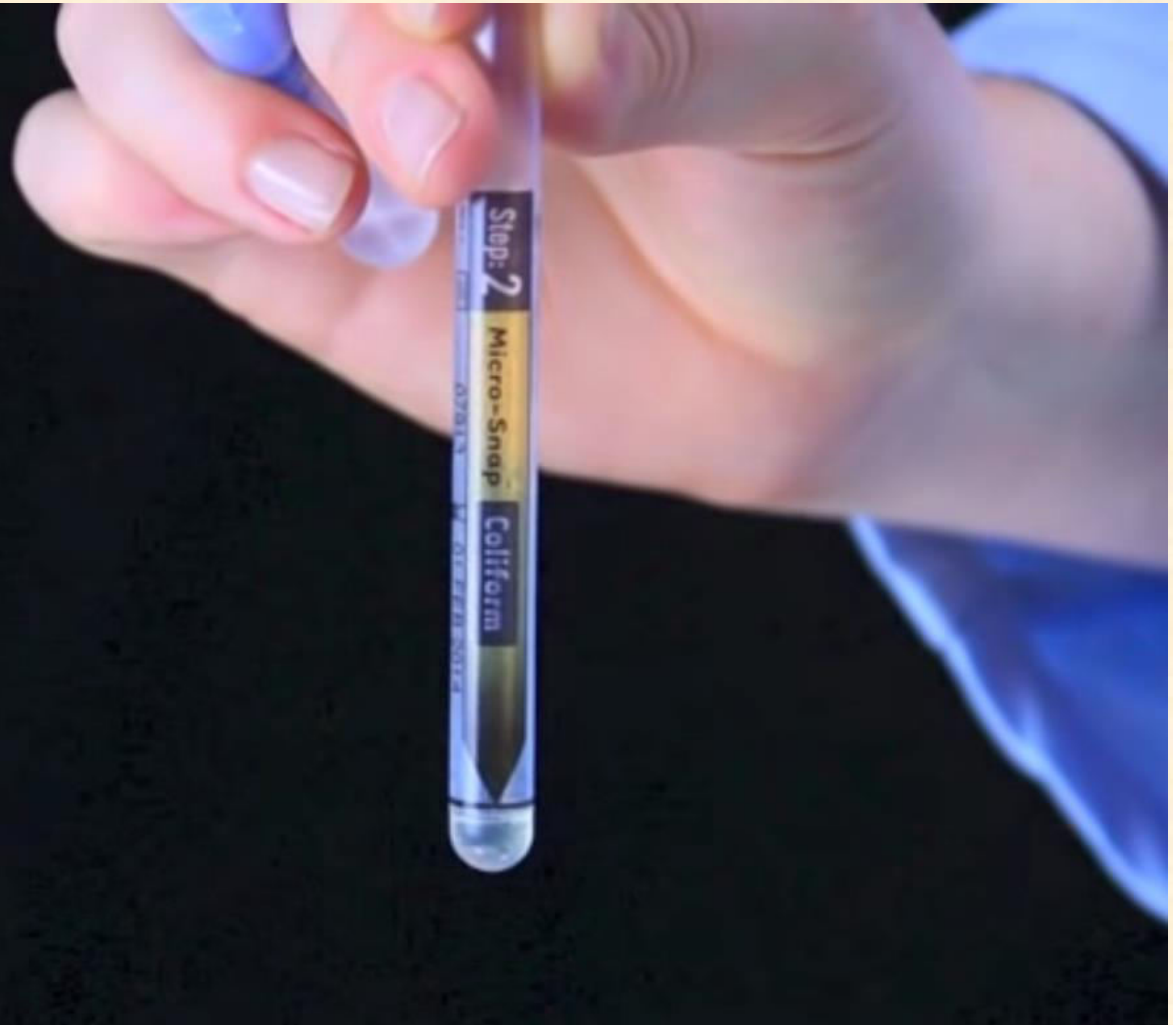Pipette enriched sample
to fill line.

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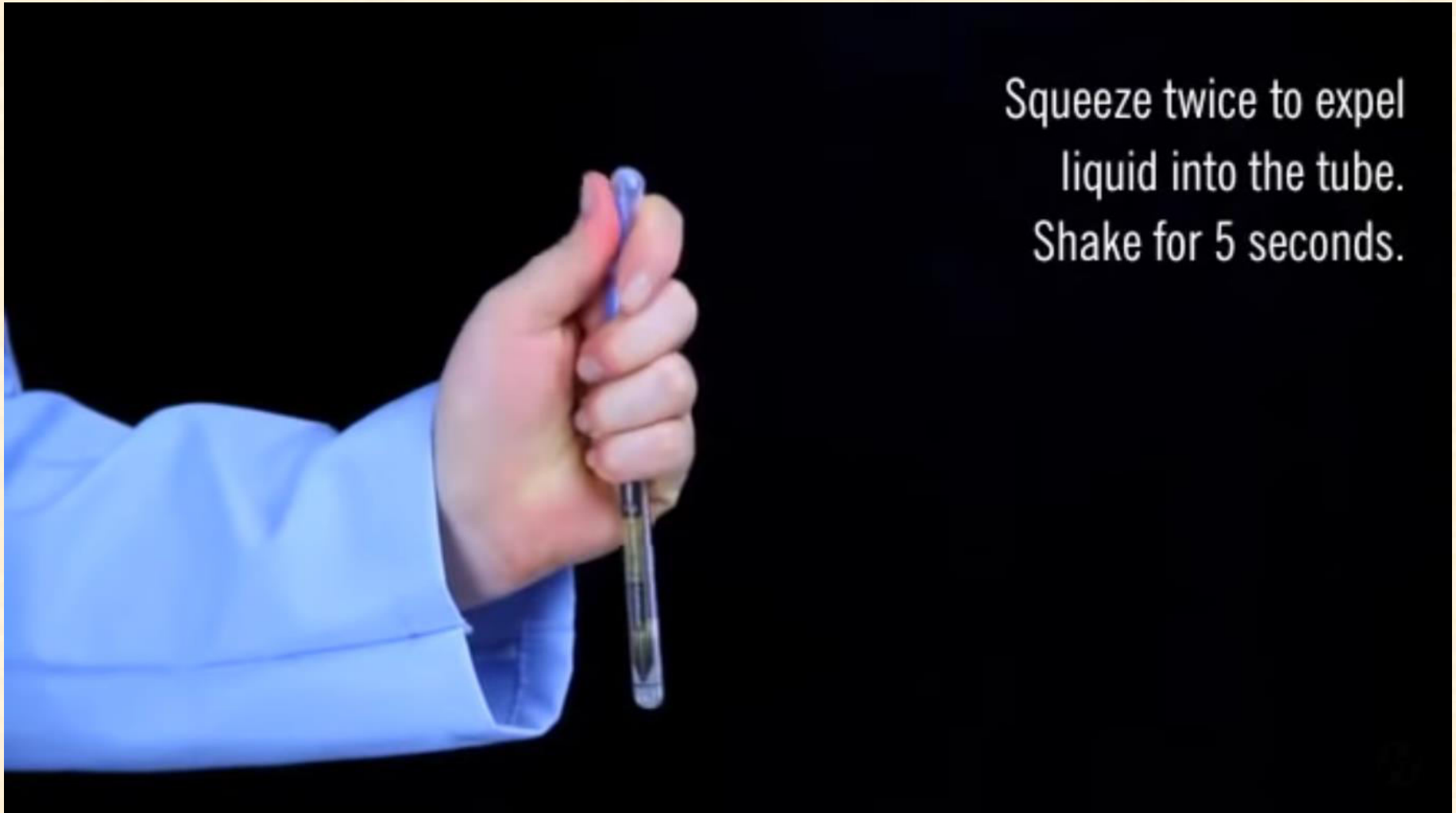
Replace the bulb on the
MicroSnap Detection Device.

Letörés


Break the Snap-Valve
by bending forward
and backward.

Letörés

Squeeze twice to expel
liquid into the tube.
Shake for 5 seconds.

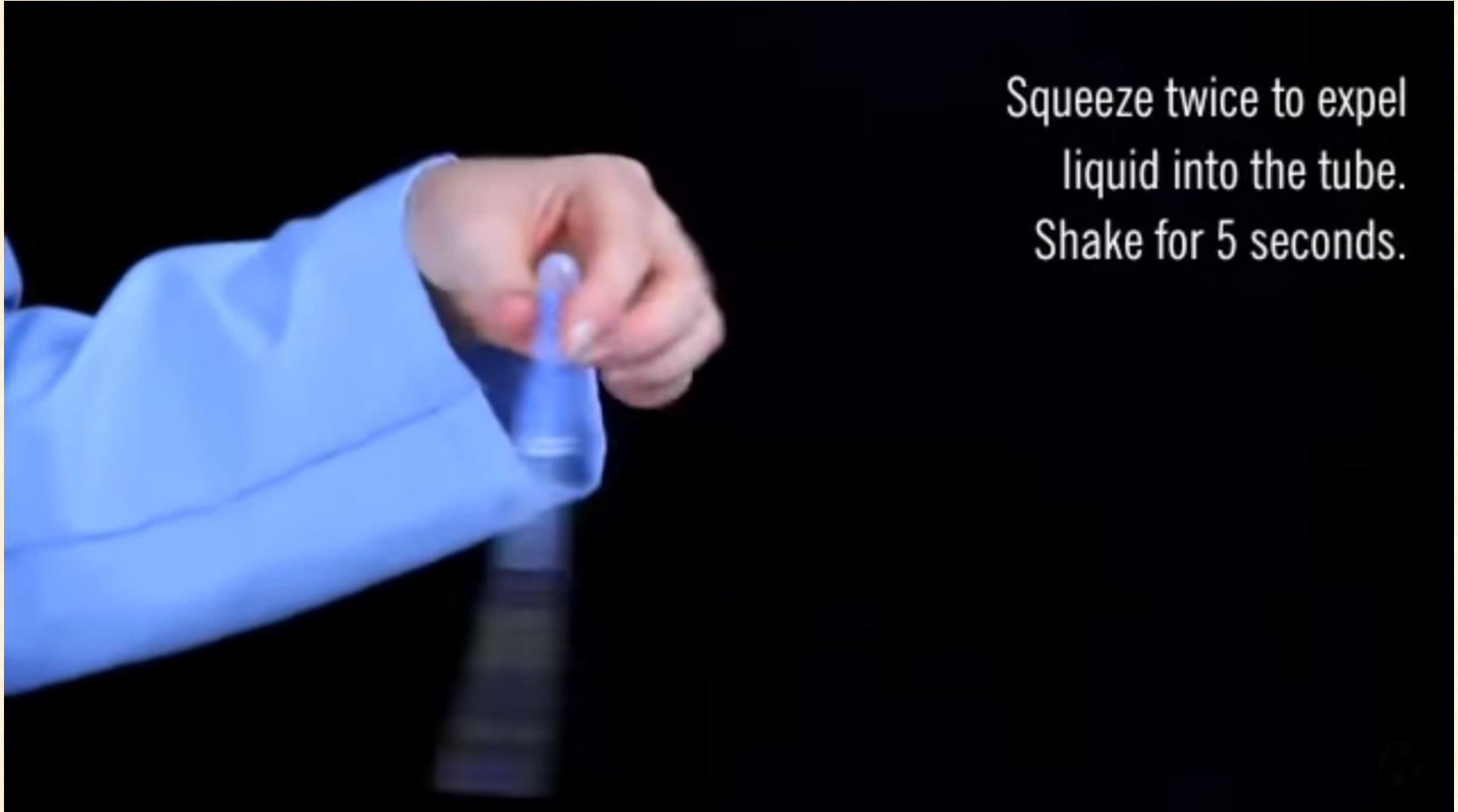Rázás

Squeeze twice to expel
liquid into the tube.
Shake for 5 seconds.

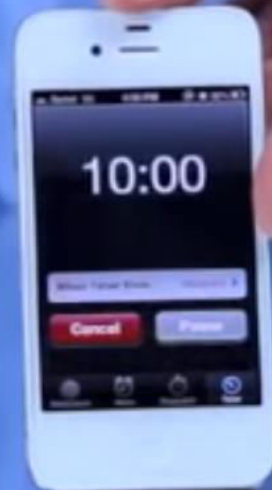
Inkubálás

Incubate **MicroSnap Detection Device**
for 10 minutes at $37^{\circ} \pm 1^{\circ}\text{C}$.

Inkubálás – 10 perc

Incubate **MicroSnap Detection Device**
for 10 minutes at $37^{\circ} \pm 1^{\circ}\text{C}$.

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ATP mérés

ATP mérés

Insert **MicroSnap** Detection Device
into **EnSURE** and initiate the
measurement.

ATP mérés

Refer to the kit insert
to interpret the results.

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
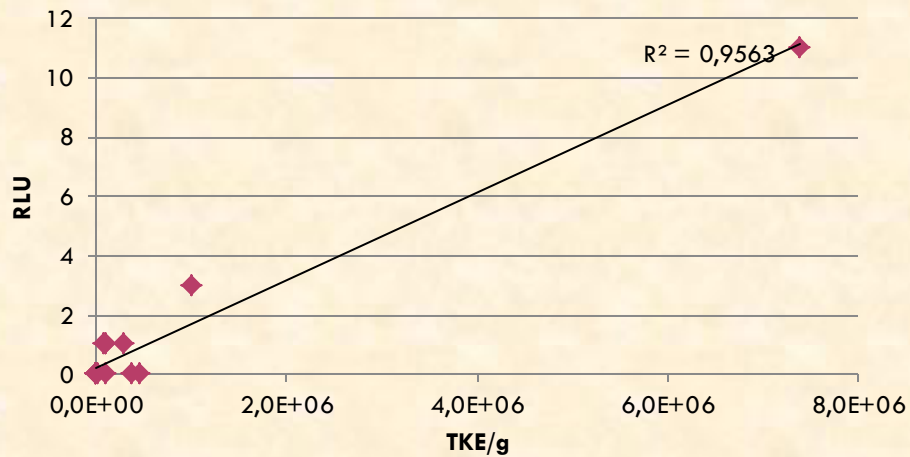
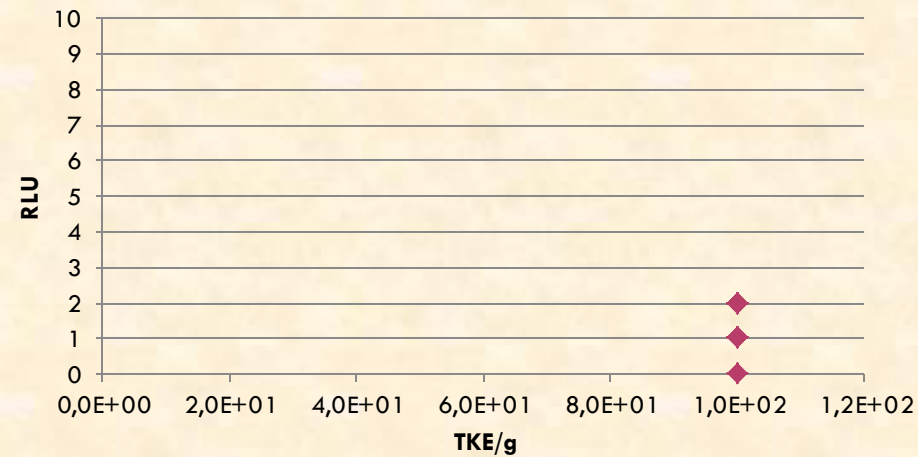
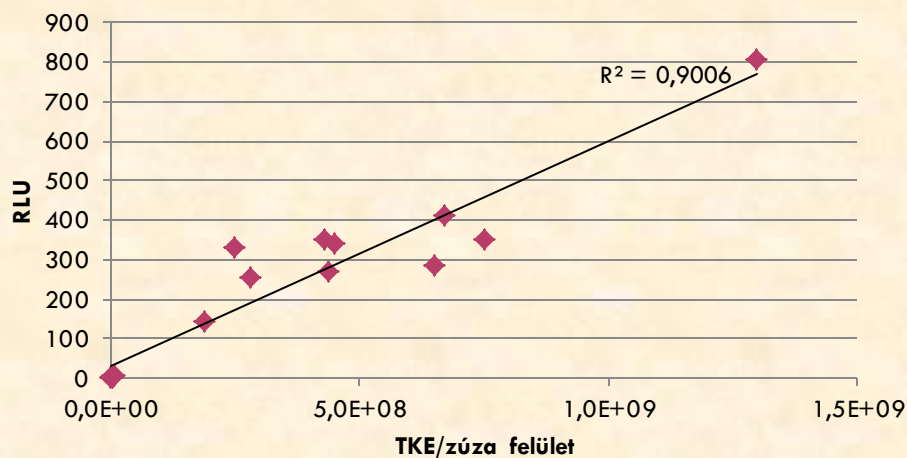
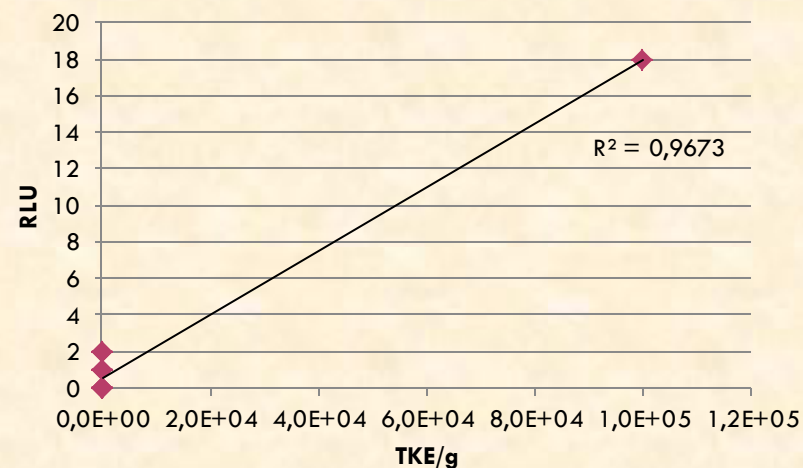
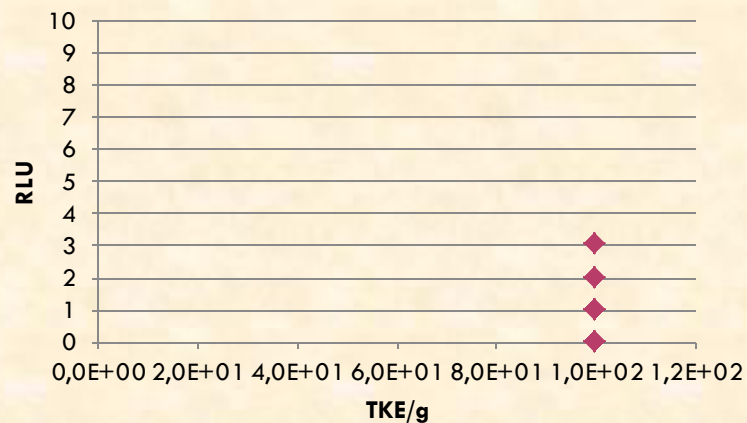
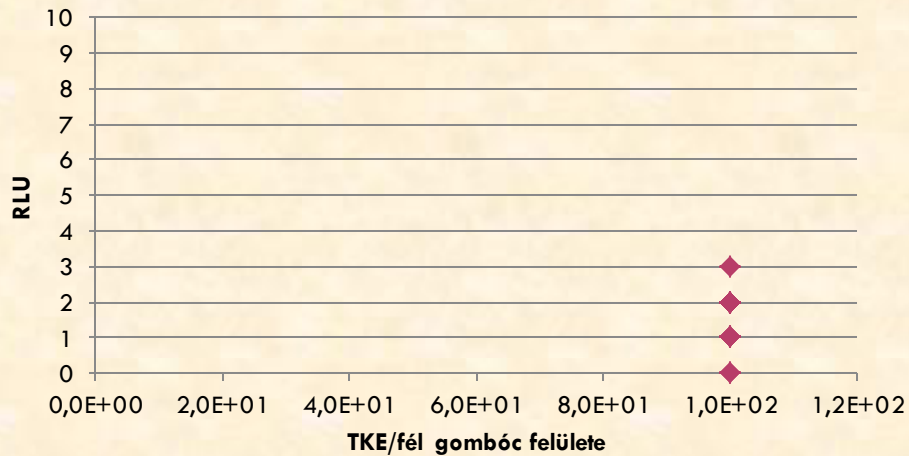
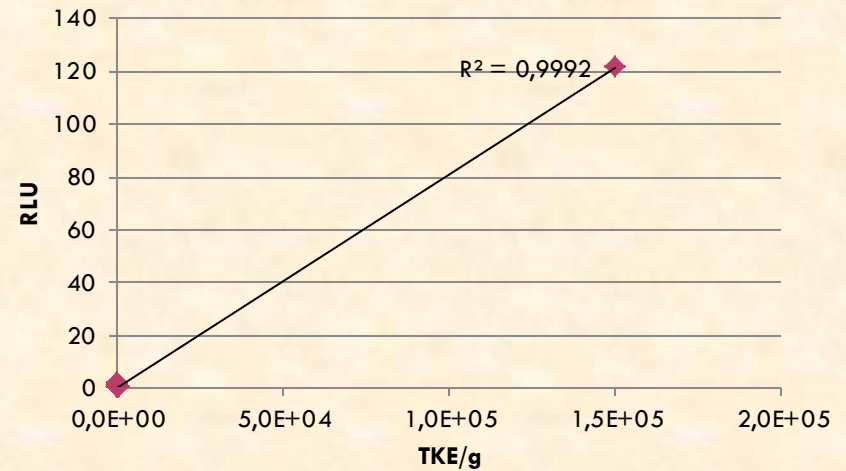
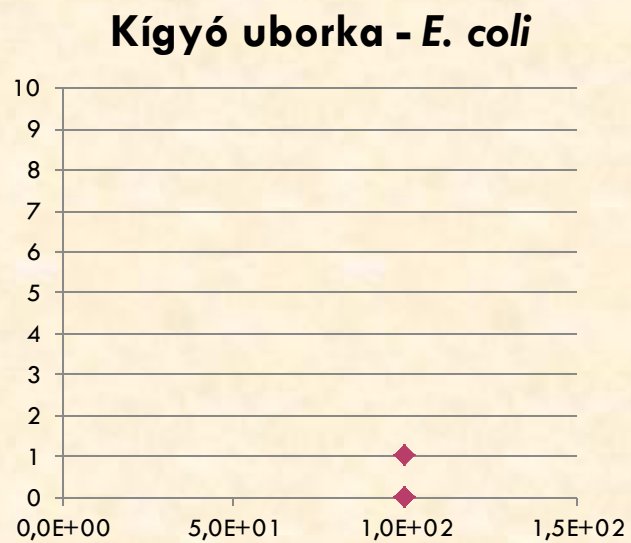
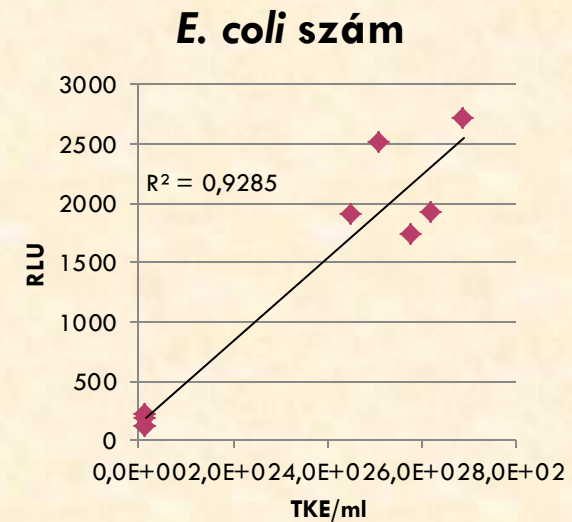
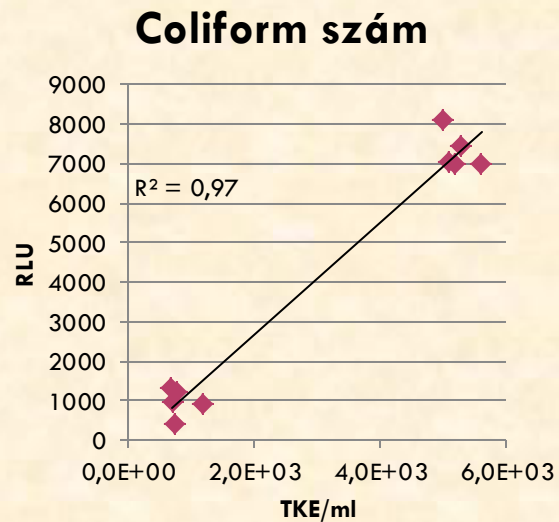
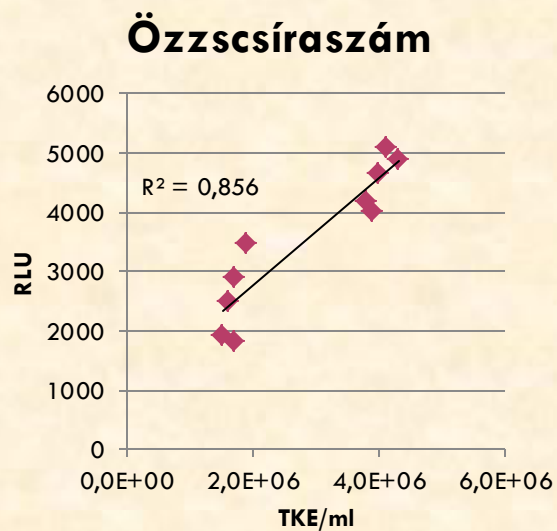
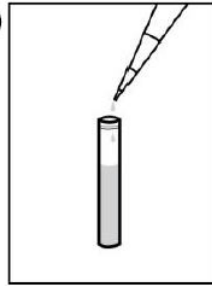
Enrichment of Milk Sample

1


1. Allow milk sample and broth to equilibrate to 22 – 26 °C. Add 1mL milk sample directly to Enhanced Nutrient Broth vial.

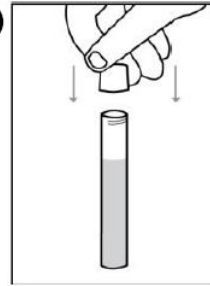2


2. Replace cap and tighten to secure.

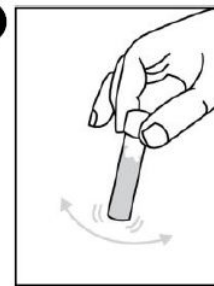3

3. Shake by hand mixing or vortexing for 10 seconds.

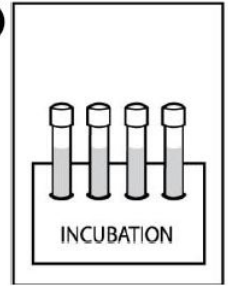4

4. Incubate at 37 ± 0.5 °C for 8 hours.

Step: 2

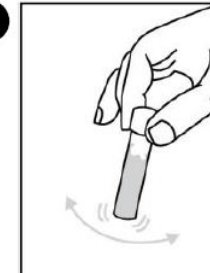
Detection / Measurement

1


1. Allow Detection Device to equilibrate to room temperature. Shake to bring liquid in tube to bottom of tube.

2


2. Shake by hand mixing or vortexing for 10 seconds.

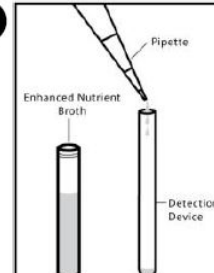3

3. Aseptically transfer 0.1mL enriched sample from Enhanced Nutrient Broth vial to Detection Device.

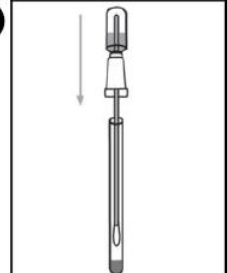4

4. Reassemble device to original state.

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Enterobacteriaceae vizsgálat

MicroSnap™ | *Enterobacteriaceae*

MicroSnap™ EB is a rapid test for detection and enumeration of *Enterobacteriaceae* (EB) bacteria. The test uses a novel bioluminogenic test reaction that generates light when enzymes that are characteristic of EB bacteria react with specialized substrates to produce light. The light generating signal is then quantified in the EnSURE™ luminometer. Depending on the required level of detection, MicroSnap EB is able to give results in the same day.

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